

PheS YEG plates

YEG medium: 0.5% yeast extract

1% NaCl

0.4% glucose

supplement with 16 mM p-Cl-Phe for YEG-pcp.

pcp is not soluble before you autoclave.

***** DO NOT AUTOCLAVE THE GLUCOSE, ADD AFTER FROM STERILE STOCK****

YEG-agar: 5 g yeast extract

10 g NaCl

2 g D,L-p-Cl-Phe

15 g agar

1 L H₂O

AUTOCLAVE

add 0.4% glucose, antibiotics

http://openwetware.org/wiki/Sauer:Making_yeg_plates

rpsL Plates:

LB medium + 15g agar.

AUTOCLAVE

Add 100µg/mL streptomycin.